

BUFFER RECIPES:

10X PBS/Glycine:

38.0 g NaCl
9.38 g Na₂HPO₄
2.07 g NaH₂PO₄
37.5 g Glycine

10X IF Wash:

38.0 g NaCl
9.38 g Na₂HPO₄
2.07 g NaH₂PO₄
2.5 g NaN₃
5.0 g BSA
10.0 ml Triton-X 100
2.05 ml Tween-20

Bring both solutions up to a total volume of 500 ml, pH 7.4 . Dilute to 1X solutions before use.